

Measurements with PRIMELAB

- 1) PRESS START BUTTON
- 2) RIGHT BUTTON: TEST
- 3) READ DESCRIPTION (OR SKIP)
- 4) CHOOSE ACCOUNTS: Default (or other), PRESS OK
- 5) CHOOSE TEST: LEGIONELLA, PRESS OK
- 6) 147-LEGIONELLA, PRESS OK
- 7) With the arrow choose ONLY ZERO AND TEST, PRESS OK
- 8) Introduce the VIAL (CONTROL), PRESS ZERO (WITH LIGHT PROTECTION)
- 9) Introduce the VIAL (SAMPLE), PRESS TEST (WITH LIGHT PROTECTION)
- 10) READ RESULTS